EN**Adding additional licenses with Windows Vista or newer****General**

Since Windows Vista, Microsoft has introduced the UAC (User Account Control), which permits running tasks either as non-administrator or as administrator (without changes of user, switching off or similar).

<http://technet.microsoft.com/en-us/library/cc709691%28WS.10%29.aspx>

<http://technet.microsoft.com/en-us/magazine/2007.06.uac.aspx>

This function can however cause difficulties with client/server installations for the programs **tiamo**, **tiBase** and **MagIC Net**.

Problem description

One would like to add additional licenses for the program; the license code for an additional 5 licenses, for example, is added (on the server) in the program part **Configuration** in the dialog window **Program administration** (accessible via menu item **Tools ▶ Program administration**, tab **Licenses**).

The file **license.mlic**, which contains the standard license code and the additional license codes, is stored in the bin directory of the program installation (**C:\Program Files\Metrohm\Program name\bin**).

If one adds the 5 licenses and then looks at this file **license.mlic**, only the standard license code will be listed (for 3 licenses).

If one highlights the file **license.mlic** and presses **Compatibility files** in the bar, then one will arrive in the directory, **C:\Users\User\AppData\Local\VirtualStore\Program Files\Metrohm\Program name\bin** ('User' being the abbreviation (UID) for the user logged in on the operating system). The file **license.mlic** with the license codes for the standard licenses and the 5 newly added licenses is to be found here, i.e. the additional licenses are stored in the **VirtualStore** of the Windows user who was logged in on the PC at the time the additional licenses were added, only.

On the server, the program recognizes all of the licenses; the operating system adds the file **license.mlic** from the directory **C:\Users\User\AppData\Local\VirtualStore\Program Files\Program name\bin** to the detail list of the directory **C:\Program Files\Metrohm\Program name\bin**.

For the Client/Server functionality of the database server, however, the file **license.mlic** in the directory **C:\Program Files\Metrohm\Program name\bin** is relevant, and here only the basic license is listed. Therefore problems are to be anticipated when adding additional clients (more than the 3 clients present in the default settings).

Remedy**▪ For new installations**

To enter additional licenses, the program must be started as administrator (position the mouse on the Program icon press down the right-hand mouse button – select **Run as administrator** there). If one then adds the license code for additional licenses in the usual way, then the file **license.mlic** in the directory **C:\Program Files\Metrohm\Program name\bin** will contain both license codes.

▪ For existing installations

First of all, the file **license.mlic** in the directory **C:\Users\User\AppData\Local\VirtualStore\Program Files\Metrohm\Program name\bin** must be deleted. This file must not be present in the **VirtualStore** of any user – a check must be made to make sure of this, and any such files which may be found must be deleted. Afterwards, the program must be started as administrator (position the mouse on the Program icon and click the right mouse button – select **Run as administrator**). If one then adds the license code for additional licenses in the usual way, then the file **license.mlic** in the directory **C:\Program Files\Metrohm\Program name\bin** will contain both license codes.

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