

1 Purpose

This document describes the fixed bugs and problems as well as improvements in the following new software versions in comparison with the previous versions.

- Vision Air
New: 2.0.5
Old: 2.0.4
 - Vision Air Routine non-pharma and pharma
New: 2.0.5.5
Old: 2.0.4.11
 - Vision Air Manager Local non-pharma and pharma
New: 2.0.5.3
Old: 2.0.4.6
 - Vision Air Manager Network non-pharma and pharma
New: 2.0.5.2
Old: 2.0.4.8

2 New features

Vision Air

Topic	Description
Instrument calibration	Instrument calibration of the DS2500 Solid Analyzer in Vision Air.

Vision Air Pharma

Topic	Description
Instrument calibration	Instrument calibration of the DS2500 Solid Analyzer in Vision Air.

3 Improvements

None

4 Fixed bugs

Vision Air

Topic	Description
Updated reports - Wavelength certification reports	In some cases, the wavelength certification report could not be created. This has been fixed.
Updated reports - Sample details report	Only 1 qualification result is displayed if hierarchical operating procedures with multiple products are used. In previous versions the hierarchical structure was not followed. Qualification results for all products were displayed. This has been fixed.

Vision Air Pharma

Topic	Description
Updated reports - Wavelength certification reports	In some cases, the wavelength certification report could not be created. This has been fixed.

5 Known bugs

Vision Air

Topic	Description
Wavelength certification test	The wavelength certification test does not use the correction files from the USB flash drive. Instead, Vision must be used to perform the wavelength certification test.

Vision Air Pharma

Topic	Description
Wavelength certification test	The wavelength certification test does not use the correction files from the USB flash drive. Instead, Vision must be used to perform the wavelength certification test.
Signing rights	The user roles System Administrator and Instrument Manager in Vision Air Local Pharma share the same signing rights. Activating signing rights for one of these user roles activates the signing rights automatically for both user roles.

6 Compliance

The current software version does not contain any modifications that affect conformity of Vision Air regarding 21 CFR Part 11 and GAMP.