

1 Purpose

This document describes new functions, the changes, and the improvements in the following firmware version:

- 57.1008.0018 for OS version 007

In comparison with the firmware version:

- 57.1008.0017 for OS version 007

2 New functions

The following new functions have been integrated:

- Initial adjustment for servo motor at time of device production
- Detection of main board version in relation to initial adjustment

3 Improvements

The following improvements were made:

- Optimizations for tests in production
- After reset, the statistics start with 1 instead of with 0
- Various improvements

4 Fixed bugs

The following bugs were fixed:

- The endpoint of the titration is reset at the time of a recalculation (TITR-20130).
- Problem with equivalence point calculation (TITR-21803).
- Eco Titrator accepts only 22 characters from Sartorius balance (TITR-23491).
- Problems with testing method C1 (SERV-25780).
- The .csv file is not saved to the USB flash drive (Defect 105131).
- Problems with "Determination time" in the user-defined pressure (Defect 109741).
- Error in decimal places with user-defined pressure (Defect 109743).

